

Residual highly polar pesticides in beer

Jan Stenzler

Application specialist LCMS

Product group LCMS

Shimadzu Deutschland GmbH

ISEAC39, July 19-22, 2016 - Hamburg, Germany

Glyphosate: background

Feb 2016:

- Glyphosate residues found in 14 german beers significant above the maximum residue limit of 0,1 µg/l for german drinking water
- Method: ELISA (LOD in beer 0,075 µg/l)
- ADI: 0,3 mg/kg body weight
- Federal Institute for risk assessment (BfR):
no risk < 1000 l/day





Glyphosate: exemplary methods

- Vegetables: EURL-SRM Quick Polar Pesticides Method (QuPPE)
- Mother's milk: BfR-method for Glyphosate (J. Agric. Food Chem. 2016, 64, 1414–1421), LOQ 1 ng/ml
- Environment: PROMOTE project (UFZ)





Beer is a difficult matrix due to

-  ion suppression for highly polar compounds
-  interferences during derivatisation with FMOC




Highly polar compounds such as glyphosate...


-  ...have a low mass and less product ions
-  ...are amphoteric




Derivatisation with FMOC

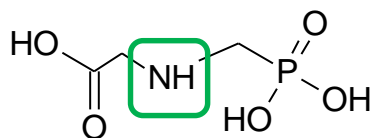


Glyphosate

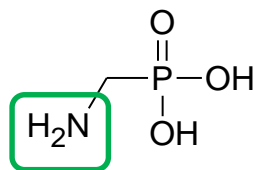

AMPA

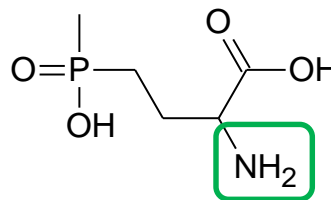


Glufosinate

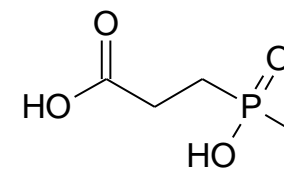

MPPA

Analyte	With derivatisation	Without derivatisation
Glyphosate	Yes	Yes
AMPA	Yes	Not possible
Glufosinate	Yes	Yes
MPPA	Not possible	Yes

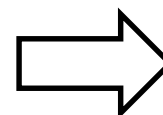

Two workflows

[1] Residual highly polar pesticides in beer: A comparative study about sample pretreatment with and without derivatization, Poster 11th EPRW, Limassol, Cyprus, May 2016

Sigrid Baumgarten¹; Anja Grüning¹; Julia Sander¹; Rebecca Kelting¹

¹ Shimadzu Europa GmbH, Duisburg, Germany

Two workflows

With derivatisation

500 μ l supernatant

+ 25 μ l EDTA-borate buffer

+ 75 μ l FMOC

60 min incubation at 50 °C

30 μ l 0,2 % phosphoric acid

+ 125 μ l water

Without derivatisation


1 ml beer + 1 ml methanol, vortex

centrifuge, 15 min, 12000 rpm

500 μ l residue

Multiplexing with Nexera MX and LCMS-8060

Nexera MX

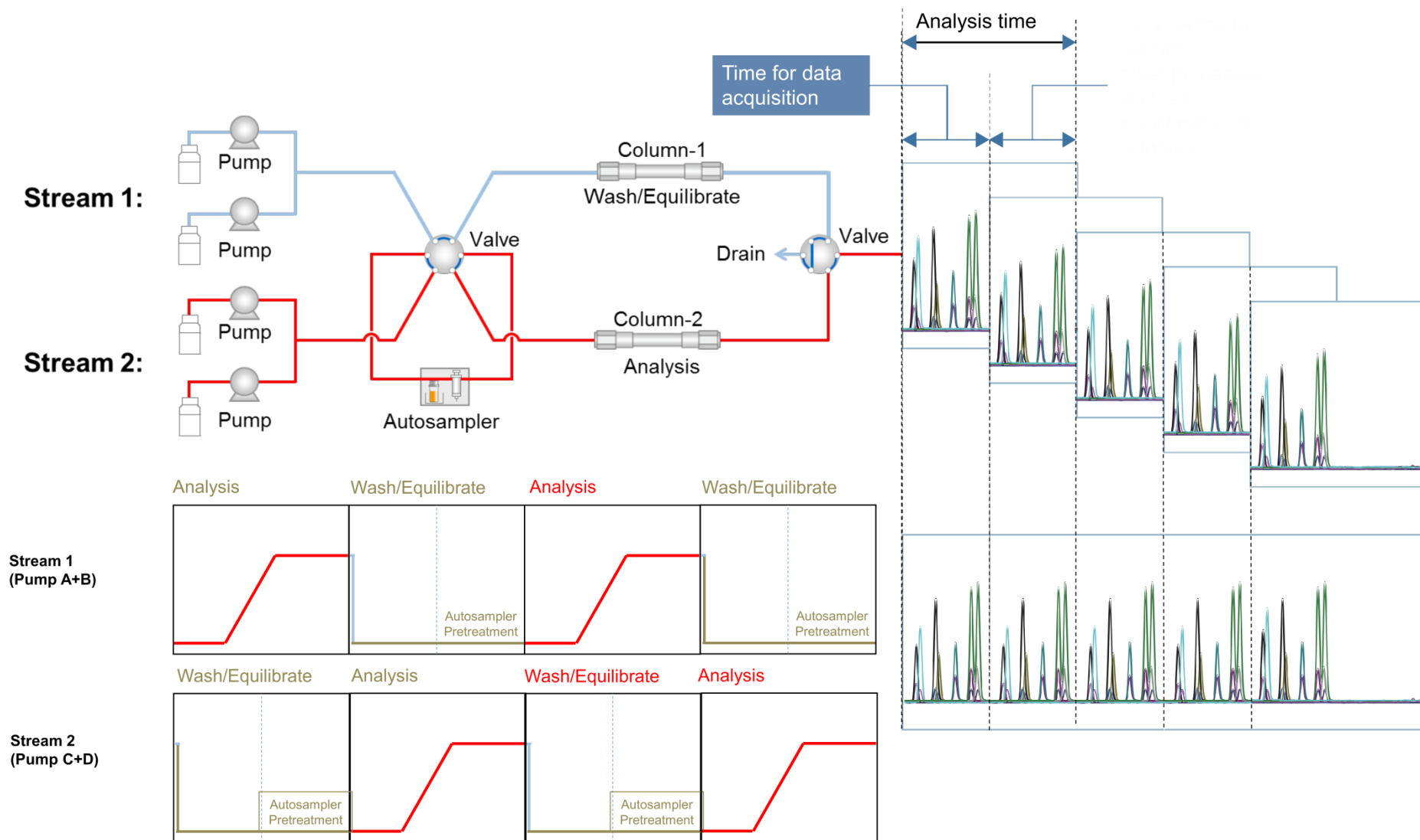

LC-MS/MS parameters

	Compared	Final
Columns, without derivatisation [2]	Obelisc R, SM-C18, SW-C18, ZIC-HILIC, ZIC-cHILIC, Hypercarb and Synergi Fusion C18	Hypercarb (100 mm x 2.1 mm, 5 µm)
Mobile phases, without derivatisation [2]	Different concentrations of ammonium acetate, ammonium formate, ammonium hydroxide, formic acid and acetic acid. Methanol, acetonitrile	1% acetic acid, methanol + 1% acetic acid
Columns, with derivatisation [1]	Raptor Biphenyl, Raptor C18 and other C18-columns	Raptor C18 (100 mm x 2.1 mm, 2.7 µm)
Mobile phases with derivatisation [3]	Shimadzu application note	5 mM ammonium acetate Acetonitrile
Ion Source Parameters, for both [1]	Drying Gas, Nebulising Gas, Heating Gas, ESI- Voltage, Desolvation Line Temperature, Heat Block Temperature, Interface Temperature	DG 3.00 l/min, NG 3.00 l/min, HG 15.0 l/min, ESI + 4 / - 3 kV, DLT 150°C, HBT 400 °C, IFT 325 °C



[2] Highly Polar Pesticide Analysis in Food Samples by LC-MS/MS, WP-351, ASMS 2014, Poster WP-351

David R. Baker¹, Mikaël Levi², Eric Capodanno³



¹Shimadzu, Manchester, UK ²Shimadzu, Marne-La-Vallée, France ³Phytocontrol, Nimes, France

[3] Application News 120, Analysis of Glufosinate, Glyphosate, and AMPA in Drinking Water Using a Triple Quadrupole LC/MS/MS System, Nov. 2015

Results: calibration and QC samples

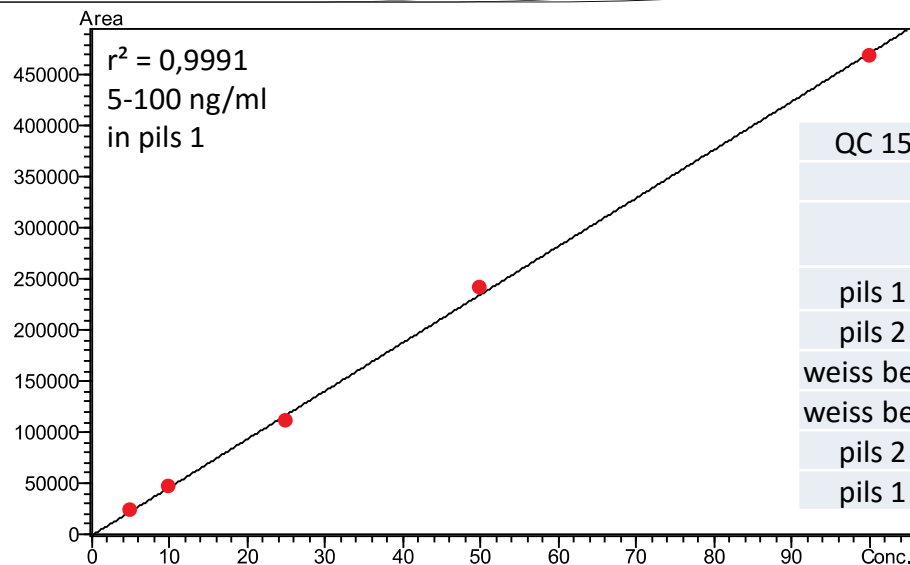
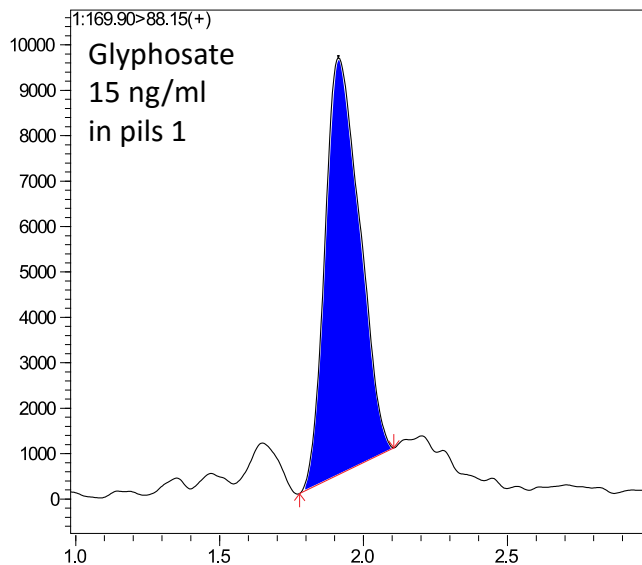


QC 15 ng/mL, spiked in beer		
Glyphosate		
	Conc.	Accuracy [%]
pils 1	15.81	105.4
pils 2	17.67	117.8
weiss beer	17.3	115.4
weiss beer	17.48	116.6
pils 2	15.62	104.1
pils 1	16.57	110.5

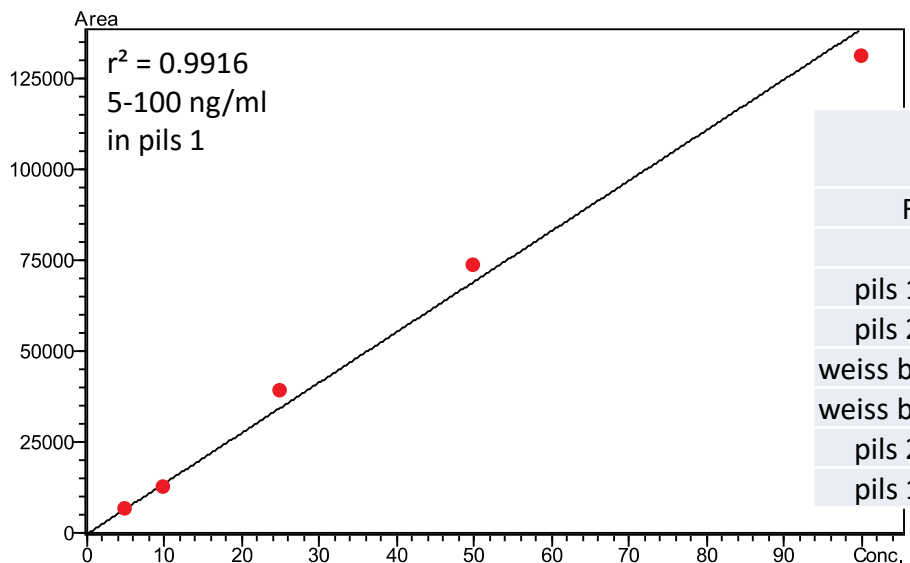
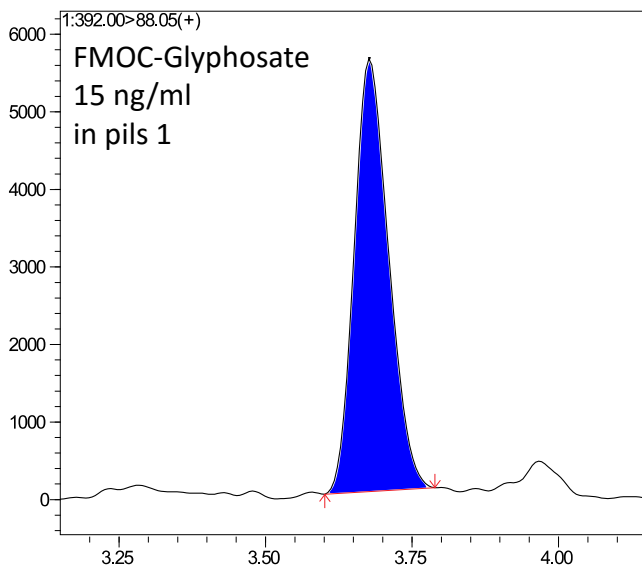


QC 15 ng/mL, spiked in beer		
FMOC-Glyphosate		
	Conc.	Accuracy [%]
pils 1	18.28	121.8
pils 2	13.04	86.9
weiss beer	16.68	111.2
weiss beer	13.32	88.8
pils 2	17.1	114
pils 1	14.42	96.1

Results: calibration and QC samples

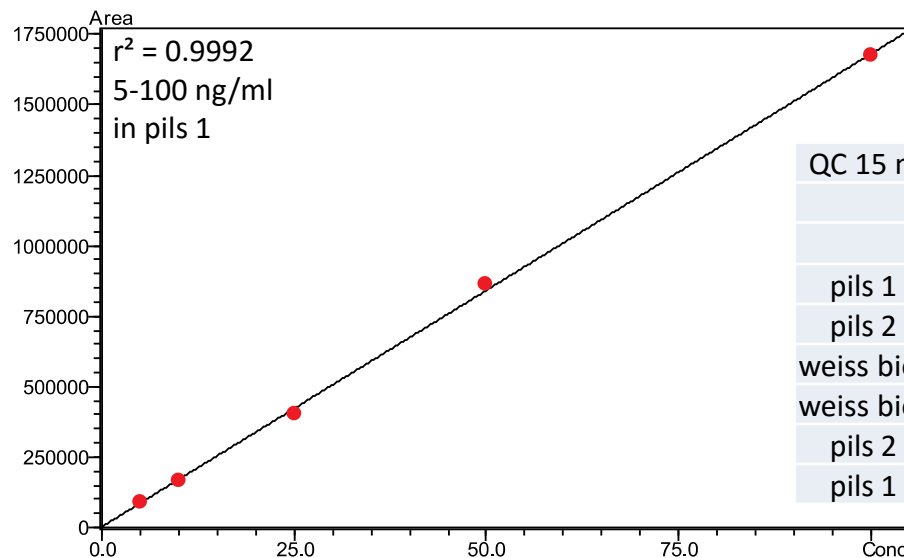
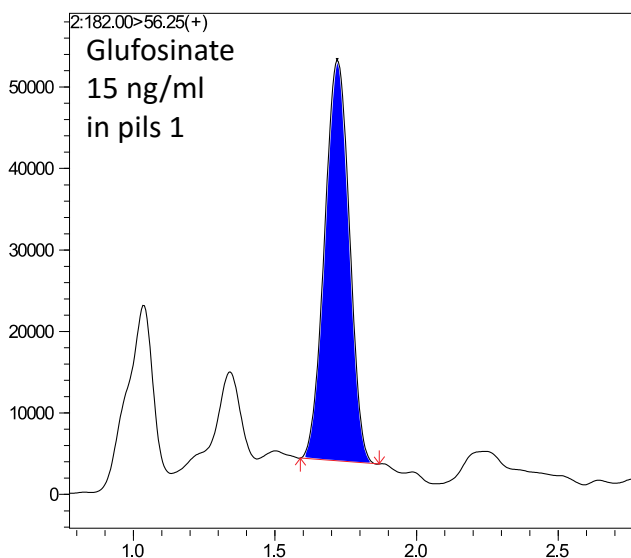


QC 15 ng/mL, spiked in beer Glufosinate		
	Conc.	Accuracy [%]
pils 1	17.12	114.1
pils 2	17.07	113.8
weiss beer	17.85	119
weiss beer	17.63	117.6
pils 2	15.35	102.3
pils 1	14.81	98.7

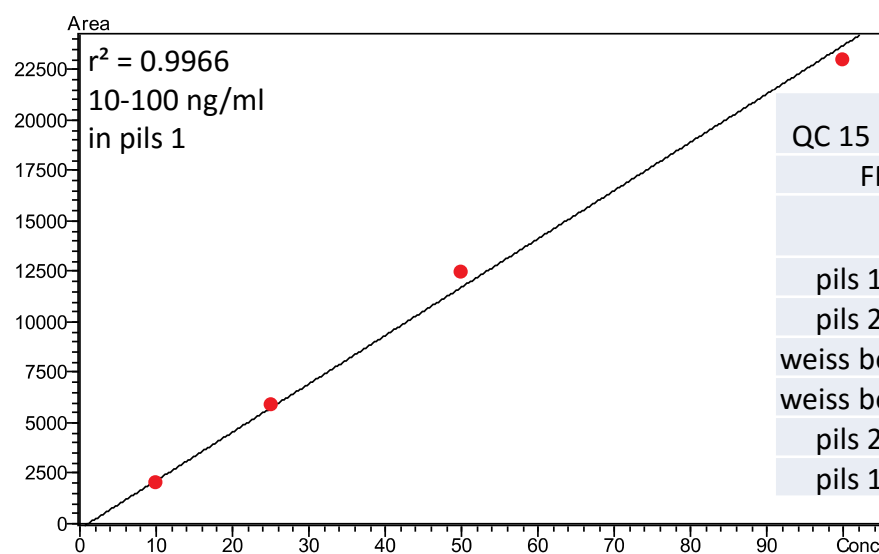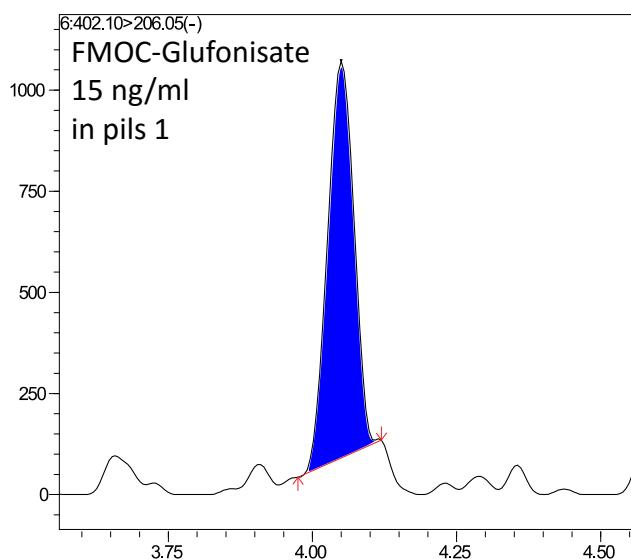



QC 15 ng/mL, spiked in beer FMOC-Glufonisate		
	Conc.	Accuracy [%]
pils 1	18.97	126.4
pils 2	15.58	103.9
weiss beer	19.77	131.8
weiss beer	16.27	108.5
pils 2	14.14	94.3
pils 1	19.81	132.1

Results: calibration and QC samples

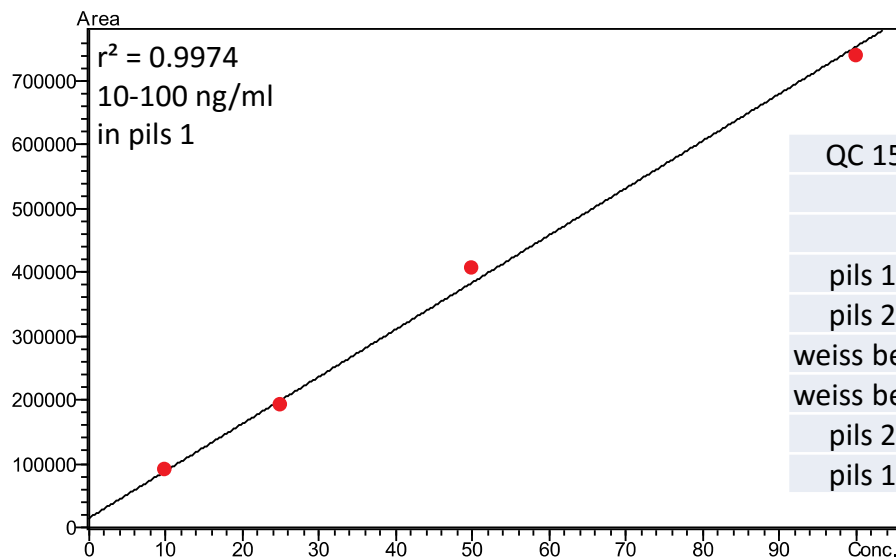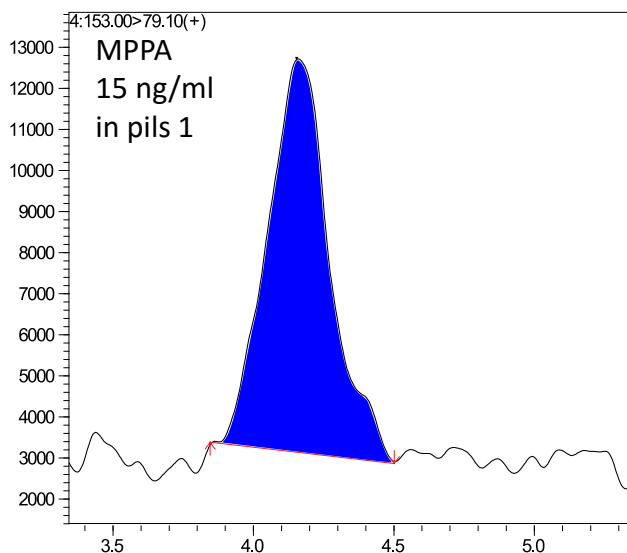

QC 15 ng/mL, spiked in beer


MPPA		
	Conc.	Accuracy [%]
pils 1	18.71	124.7
pils 2	15.23	101.5
weiss beer	17.91	119.4
weiss beer	17.09	113.9
pils 2	14.08	93.9
pils 1	18.12	120.8

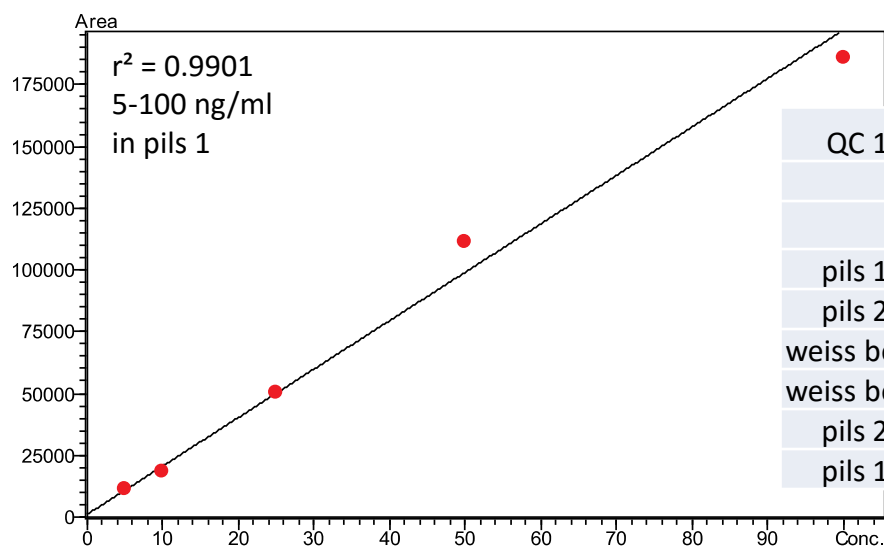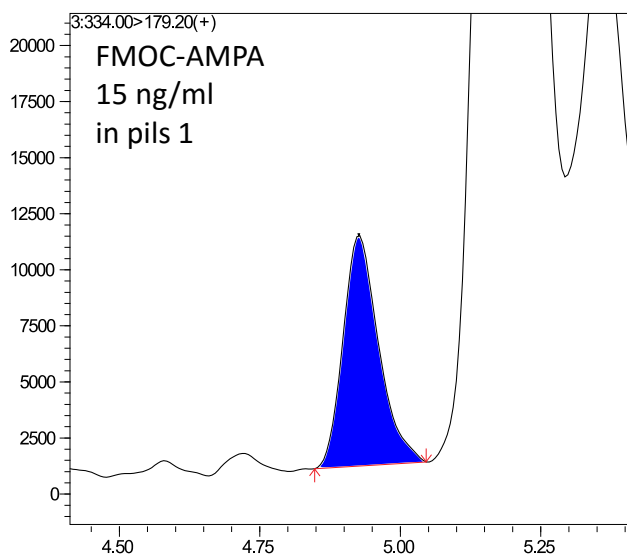

QC 15 ng/mL, spiked in beer

FMOC-AMPA		
	Conc.	Accuracy [%]
pils 1	22.7	151.4
pils 2	13.31	88.7
weiss beer	15.9	106
weiss beer	17.53	116.9
pils 2	12.85	85.6
pils 1	16.98	113.2

Results: beer samples

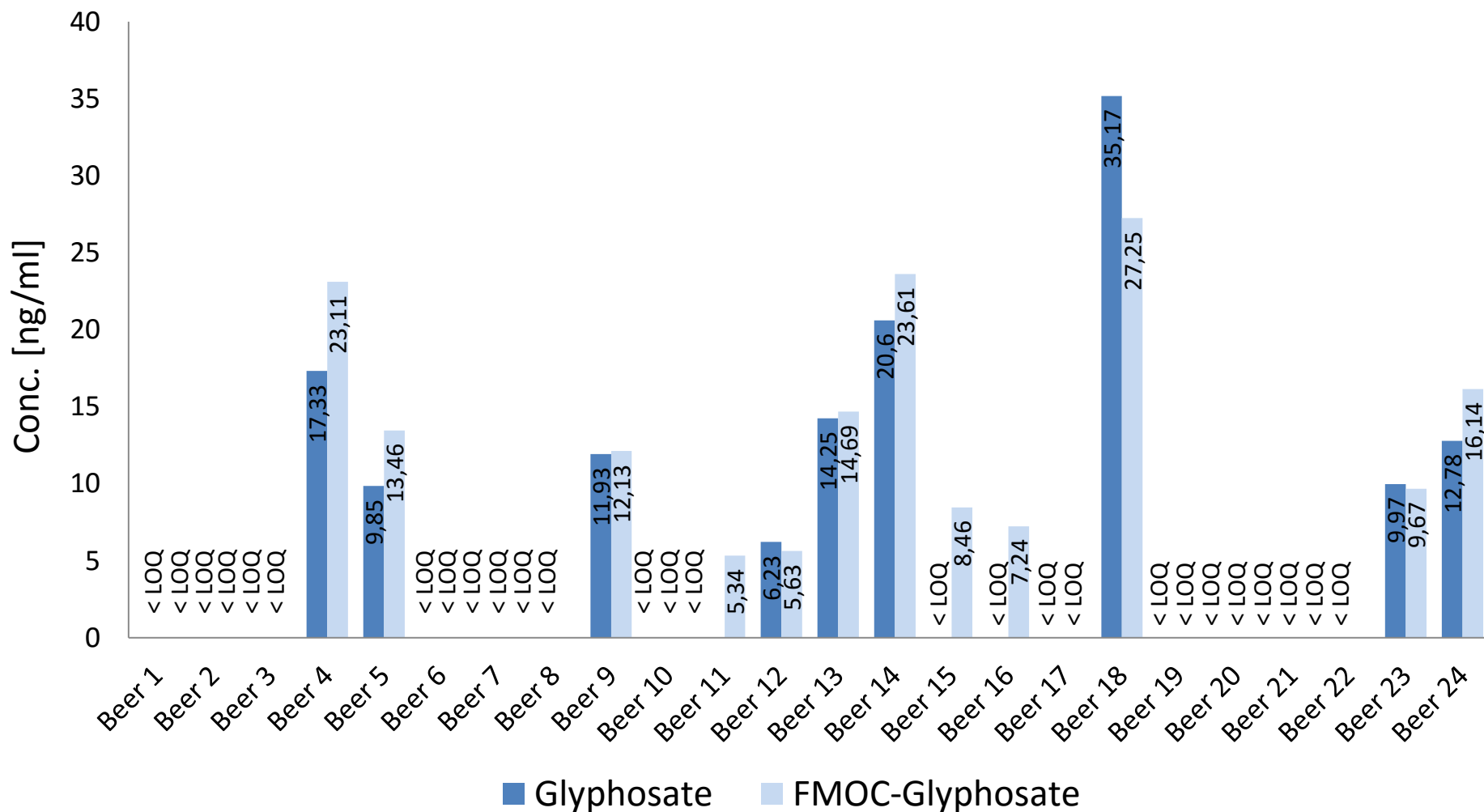

Glyphosate and AMPA in food



LCMS-8050: confirm results at low levels with fast polarity switching (5 msec)

FMOC-Glyphosate
(ESI positive)




FMOC-Glyphosate
(ESI negative)

Quantification and
identification by
standard addition

organic grain sample

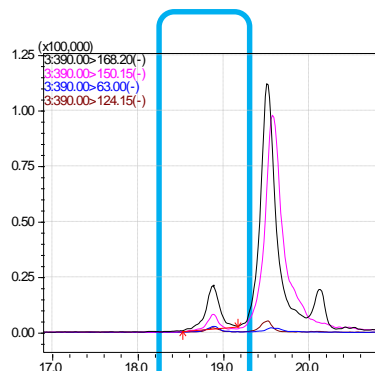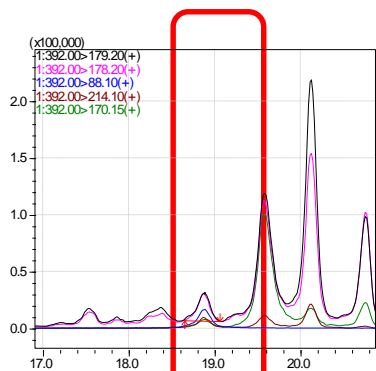

organic grain sample
+ 250 ppb

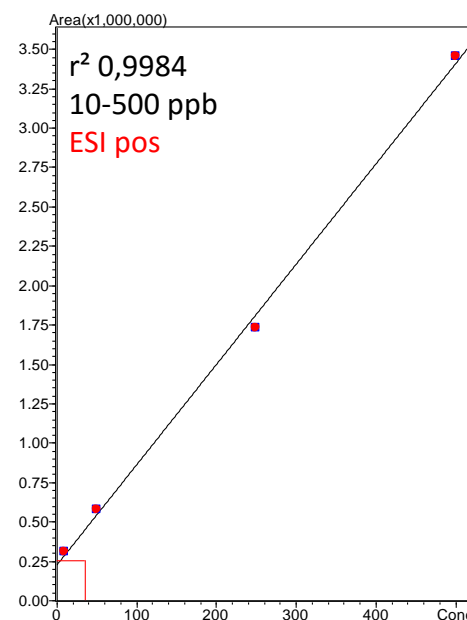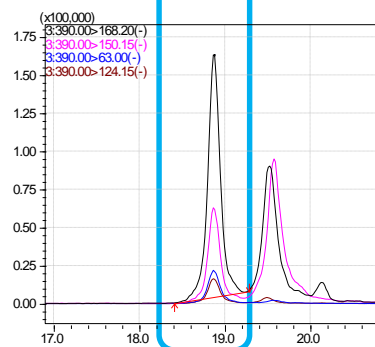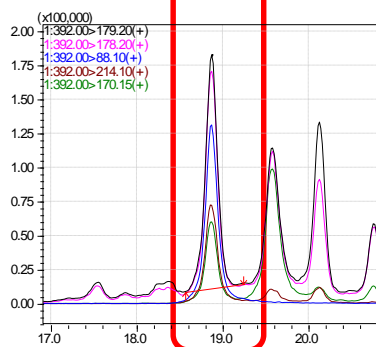

Glyphosate and AMPA in food

LCMS-8050: confirm results at low levels with fast polarity switching (5 msec)

FMOC-AMPA
(ESI positive)



FMOC-AMPA
(ESI negative)

Quantification and
identification by
standard addition

organic grain sample



organic grain sample
+ 250 ppb

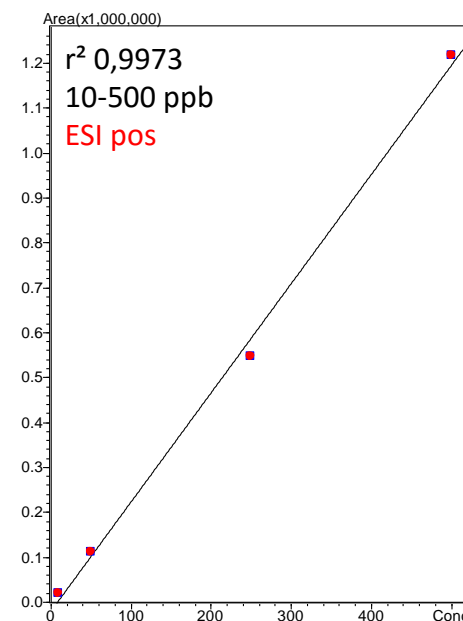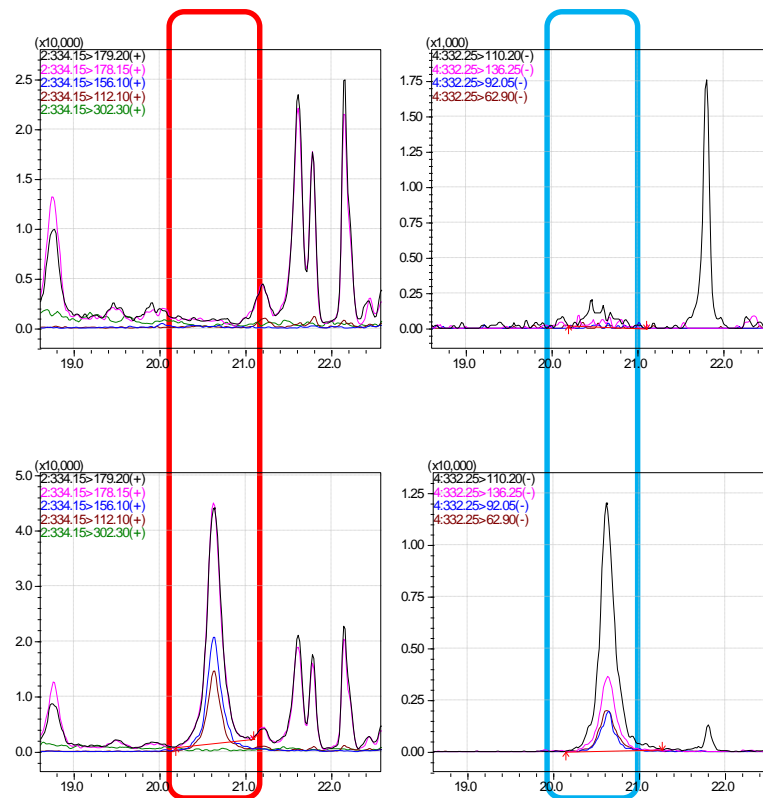

Summary

- A routine method for the determination of Glyphosate, AMPA, Glufosinolate and MPPA was developed
- LOQ: 10 ng/ml in beer
- Accurate and sensitive quantification without derivatisation
- Nexera MX allows higher sample throughput
- Application Notes available (request: info@shimadzu.eu)



Residual highly polar pesticides in beer: A comparative study about sample pretreatment with and without derivatization

Sigrid Baumgarten¹; Anja Grüning¹; Julia Sander¹; Rebecca Kelting¹

¹Shimadzu Europa GmbH, Duisburg, Germany



Highly Polar Pesticide Analysis in Food Samples by LC-MS/MS

David R. Baker¹, Mikaël Levi², Eric Capodanno³

¹Shimadzu, Manchester, UK ²Shimadzu, Marne-La-Vallée, France ³Phytocontrol, Nîmes, France

Application
News

No. **C120**

Liquid Chromatography Mass Spectrometry

Analysis of Glufosinate, Glyphosate, and AMPA in Drinking Water Using a Triple Quadrupole LC/MS/MS System

Thank you



Brand statement “Excellence in Science”

We in the Shimadzu Group have delivered products and services to enable our customers around the world to develop a diverse range of new products, to protect and improve the environment, and to improve the health and lives of mankind. This brand statement represents our sense of pride in this endeavor. It is our commitment to society and ourselves that Shimadzu remains dedicated in our pursuit of technology and accumulation of knowledge, so that we can offer even more outstanding technologies, products, and services, so as to be recognized for excellence in the field of science.